# Levenhuk 2L NG/3L NG/D2L NG

EN User Manual

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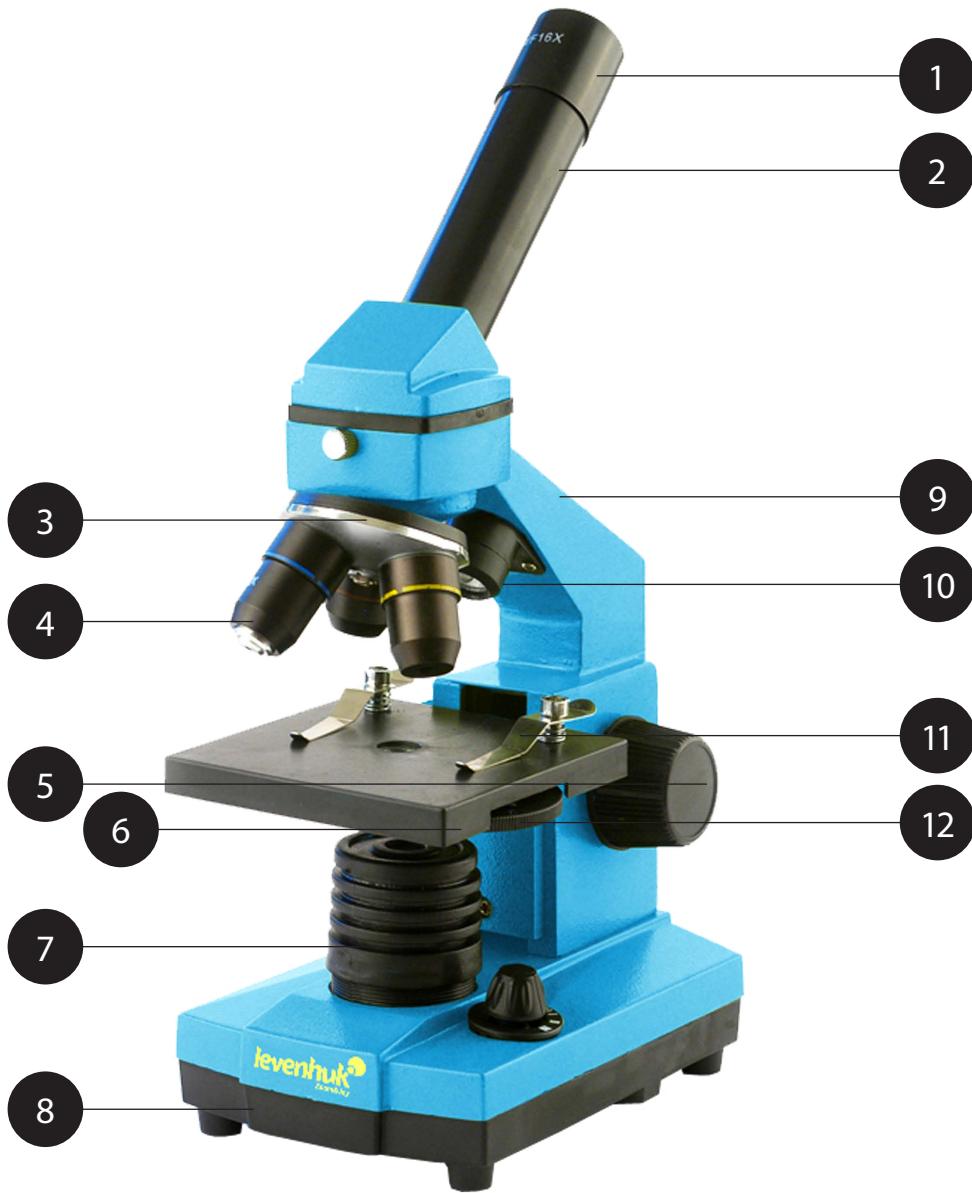
16

## General Information

The microscope is safe for health, life and property of the consumer and the environment when properly used; the microscope meets the requirements of international standards. The microscope is designed for observing transparent objects in the transmitted and reflected light using the bright field method, for biological use and teaching demonstrations.

Levenhuk D2L digital camera was exclusively designed for use with this microscope. It allows transfer of the precise image of the observed object to your PC display. Levenhuk ToupView software is included in the package; it allows viewing and editing of the received images.

## Specification

Magnification	64x–640x
Magnification of objective	4x, 10x, 40x
Eyepiece magnification	16x
Stage dimensions	90x90mm
Stage moving range	0–10mm
Power supply	110V/50 Hz AC (adapter), 3 AA batteries
Illumination (upper/lower light)	LED
Camera model	Levenhuk D2L
MAX resolution (at rest)	640x480
Megapixels	0.3
Sensor	1/4" CMOS
Pixel size	6µm x 6µm
Sensitivity, v/lux·sec @550 nm	2.0
Speed (depends on the PC model)	30 fps
Video recording	Yes
Active range	60dB
Location	ocular tube (not eyepiece)
Image format	*.bmp, *.jpg, *.jpeg, *.png, *.tif, *.tiff, *.gif, *.psd, *.ico, *.emf, etc.
Spectral range	400–650nm
Programmable options	Image size, brightness, white balance, exposure control, etc.
Slot	USB 2.0, 480 Mb/s
System requirements	Windows XP/Vista/7/8 (32-bit and 64-bit), minimum Intel Core 2 2.8 GHz, USB 2.0 port
Software	USB 2.0 driver, Levenhuk ToupView
Power supply	via USB cable
Temperature range	-30...70 °C, -86...158 °F

The manufacturer reserves the right to make changes to the product range and specifications without prior notice.

## Legend

- ① Eyepiece
- ② Monocular nosepiece
- ③ Revolving head
- ④ Objectives
- ⑤ Focusing knob
- ⑥ Stage
- ⑦ Lower illumination
- ⑧ Base stand
- ⑨ Stand
- ⑩ Upper illumination
- ⑪ Slide holders
- ⑫ Diaphragm disk
- ⑬ USB port
- ⑭ Digital camera
- ⑮ USB cable
- ⑯ Setup CD

## Package

### Levenhuk 2L NG

Microscope unit with stage, revolving head for three objectives, focusing mechanism, two illuminations	1 Pc
Monocular nosepiece	1 Pc
Objectives: 4x, 10x and 40x	1 Pc each
Eyepiece WF16x	1 Pc
AC adapter	1 Pc
AA batteries	3 Pcs
User manual and warranty slip	1 Pc each

### Levenhuk D2L NG

In addition to Levenhuk 2L NG: Levenhuk D2L digital camera:

Digital camera	1 Pc
USB cable	1 Pc
SETUP CD Levenhuk ToupView (auto setup guide, drivers, software for image visualization, fixation and processing)	1 Pc

### Levenhuk 3L NG

In addition to Levenhuk 2L NG: Levenhuk K50 experiments kit:

Prepared slides	5 Pcs
Clean slides	5 Pcs
Flasks: with pitch, sea salt, yeast, brine shrimp	1 Pc each
Brine shrimp hatchery	1 Pc
Microtome	1 Pc
Forceps	1 Pc
"Attractive microscope. Scrutinizing the microcosm" user guide	1 Pc

**Caution:** Please refer to the specifications table for the correct mains voltage and never attempt to plug a 110V device into 220V outlet and vice versa without using a converter. Remember that mains voltage in the U.S. and Canada is 110V and 220–240V in most European countries.

## Modules description and usage

Focusing mechanism located on the stand allows vertical movement of the stage. The stage is moved by focusing knob. The total focusing value is at least 10mm.

Slide holders are fixed on the stage surface. Diaphragm disk is attached to the lower part of the stage.

Revolving head holds the objectives; the objectives can be replaced by rotating the ring until it reaches the fixed position.

Objectives:

Linear magnification	Numerical aperture
4	0.1
10	0.25
40	0.65

The 40x objective is equipped with a spring casing, which protects the frontal lens from contacting the specimen during the focus adjusting.

## Using the microscope

### Unpacking and getting started

- Unpack the microscope.
- Make sure all the parts are present.
- Move the stage to the lowermost position using the focusing knob.
- Press the ON/OFF button located on the back side of the microscope base next to the power cord to turn the microscope on.

### Focusing

- Place an object on the stage and fix it with the slide holders.
- Rotating the revolving head, switch magnification to the 4x objective.
- Move the object to fix its thickest part exactly under the objective.
- Rotate the focusing knob to raise the stage until the objective is close to the specimen; keep checking the distance between the objective and the object to avoid their contact.
- Look through the eyepiece installed in the monocular head and lower the stage slowly rotating the focusing knob until you see the object image.

Such adjustment protects the frontal lens from contacting the object when you use objectives of other magnifications; though, slight refocusing might be required.

### Selecting the objective

We recommend that you start your observations with the lowest power objective; it helps to select a segment for the detailed research. Once the segment is selected, you should center its image in the microscope's field of view. This operation should be performed as precisely as possible, otherwise the required segment might fail to appear in the field of view of the higher power objectives.

## Camera

The digital camera was exclusively designed for use with this microscope.

The object image can be transferred to your PC display in its actual colors and stored in the PC memory.

Levenhuk ToupView software is included in the package; it allows viewing and editing of the received images. Supported file formats: \*.bmp, \*.jpg, \*.jpeg, \*.png, \*.tif, \*.tiff, \*.gif, \*.psd, \*.ico, \*.emf, etc.

The USB cable is used as camera power supply and for connection with PC.

Compatible OS: Windows XP/Vista/7/8 (32-bit and 64-bit), minimum Intel Core 2 2.8 GHz, USB 2.0 port.

## Care and maintenance

- Never, under any circumstances, look directly at the Sun, another bright source of light or at a laser through this device, as this may cause PERMANENT RETINAL DAMAGE and may lead to BLINDNESS.
- Take necessary precautions when using the device with children or others who have not read or who do not fully understand these instructions.
- After unpacking your microscope and before using it for the first time check for integrity and durability of every component and connection.
- Do not try to disassemble the device on your own for any reason. For repairs and cleaning of any kind, please contact your local specialized service center.
- Protect the device from sudden impact and excessive mechanical force. Do not apply excessive pressure when adjusting focus. Do not overtighten the locking screws.
- Do not touch the optical surfaces with your fingers. To clean the device exterior, use only special cleaning wipes and special optics cleaning tools from Levenhuk. Do not use any corrosive or acetone-based fluids to clean the optics.
- Abrasive particles, such as sand, should not be wiped off lenses, but instead blown off or brushed away with a soft brush.
- Do not use the device for lengthy periods of time, or leave it unattended in direct sunlight. Keep the device away from water and high humidity.

- Be careful during your observations, always replace the dust cover after you are finished with observations to protect the device from dust and stains.
- If you are not using your microscope for extended periods of time, store the objective lenses and eyepieces separately from the microscope.
- Store the device in a dry, cool place away from hazardous acids and other chemicals, away from heaters, open fire and other sources of high temperatures.
- Seek medical advice immediately if a small part or a battery is swallowed.

## Battery safety instructions

- Always purchase the correct size and grade of battery most suitable for the intended use.
- Always replace the whole set of batteries at one time; taking care not to mix old and new ones, or batteries of different types.
- Clean the battery contacts and also those of the device prior to battery installation.
- Make sure the batteries are installed correctly with regard to polarity (+ and -).
- Remove batteries from equipment that is not to be used for an extended period of time.
- Remove used batteries promptly.
- Never attempt to recharge primary batteries as this may cause leakage, fire, or explosion.
- Never short-circuit batteries as this may lead to high temperatures, leakage, or explosion.
- Never heat batteries in order to revive them.
- Remember to switch off devices after use.
- Keep batteries out of the reach of children, to avoid risk of ingestion, suffocation, or poisoning.
- Utilize used batteries as prescribed by your country laws.

## Levenhuk International Lifetime Warranty

All Levenhuk telescopes, microscopes, binoculars and other optical products, except for accessories, carry a **lifetime warranty** against defects in materials and workmanship. **Lifetime warranty** is a guarantee on the lifetime of the product on the market. All Levenhuk accessories are warranted to be free of defects in materials and workmanship for **six months** from date of retail purchase. Levenhuk will repair or replace such product or part thereof which, upon inspection by Levenhuk, is found to be defective in materials or workmanship. As a condition to the obligation of Levenhuk to repair or replace such product, the product must be returned to Levenhuk together with proof of purchase satisfactory to Levenhuk.

This warranty does not cover consumable parts, such as bulbs (electrical, LED, halogen, energy-saving and other types of lamps), batteries (rechargeable and non-rechargeable), electrical consumables etc.

For further details, please visit our web site: <http://www.levenhuk.com/warranty>

If warranty problems arise, or if you need assistance in using your product, contact the local Levenhuk branch:

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Levenhuk is a registered trademark of Levenhuk, Inc.
© 2006–2015 Levenhuk, Inc. All rights reserved.
Levenhuk, Inc. 1935 Brandon Court, Suite A-1