# Levenhuk 2ST, Levenhuk 3ST Stereo Microscopes

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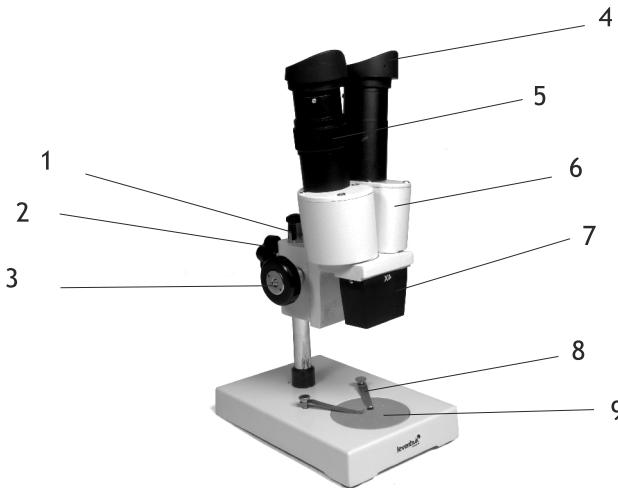
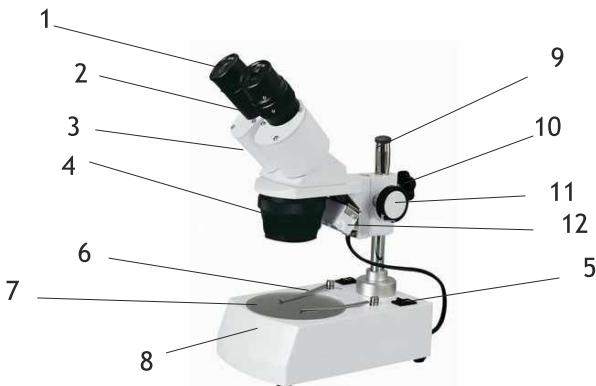
Please read this manual carefully before using your microscope for the first time.

Levenhuk 2ST and Levenhuk 3ST microscopes were designed to be used during assembly of intricate equipment, including electronics; for school experiments and scientific research. Both models may be found in schools and research institutes, at home and in manufacturing plants.

## Specifications

- Vertical binocular head (Levenhuk 2ST) or a binocular head inclined at a 45° angle (Levenhuk 3ST)
- Diopter adjustment range: +/-5D
- Interpupillary distance adjustment range: 60 mm
- Magnification: 40x (Levenhuk 2ST), 20-40x (Levenhuk 3ST)
- Objective lenses: 4x (Levenhuk 2ST), 2x, 4x (Levenhuk 3ST)
- Power source: 220V/50Hz or 110V/60Hz (optional)
- Illumination sources: Upper and lower illumination sources, 12V/10W halogen lamps (for observations in reflected or transmitted light) (Levenhuk 3ST). Natural illumination is used in Levenhuk 2ST model.

## Additional accessories

- Dark field condenser (black and white plate)
- Levenhuk 3ST: Fluorescent lamp (may be held in place with three locking screws on the side of the nosepiece). A regular lamp may be used as an alternative, for softer lighting and more comfortable observations.
- Rubber eyecups

## Using the microscope

### Environment

These microscopes perform best in dry and clean rooms - even a smallest speck of dust on the optics may reduce the quality of your views - with temperatures ranging from -5 to +40 °C (23-104 °F).

### Illumination (Levenhuk 3ST)

Plug your microscope into a power source, turn it on and choose the required illumination source. If your microscope is fitted with brightness control, you can also choose the brightness level of illumination.

### Stage inset glass

- Frosted glass stage inset may be used for observations of transparent specimens or thin slices of objects. Place it on the stage and secure it in place with the locking screw. Lower illumination source should be used during observations with this glass stage inset.
- The kit of your new microscope may also include a black and white glass stage inset. Normally, you would want to use the white side; however, if a specimen is very bright, or colored white, you should use the black side for additional contrast. Upper illumination source should be used during observations with this glass stage inset. Before installing the black and white glass make sure to remove the frosted glass from the stage.

## **Preparations**

Place a specimen in the center of a blank slide, put the slide on the stage and lock it in place with stage clips.

## **Eyecups**

For additional comfort during observations you can use rubber eyecups on both eyepieces.

## **Focusing, diopter adjustment, interpupillary distance**

Place a slide on the stage and lock it in place with stage clips. Loosen the locking knob on the microscope pole and slide the head up or down into a desired position and retighten the locking knob. Look through the right eyepiece and turn the objective lens handle until you see the specimen in your field of view. Afterwards, turn the focus knob until the image is sharp and clear. Look through the left eyepiece and rotate the diopter adjustment ring until the image is sharp and clear. Do not adjust focus during diopter adjustment! Upon completing this process, look through both eyepieces, hold them with your hands and bring them closer or move them apart until a double image turns into a single image.

## **Changing objective lenses**

Levenhuk 3ST has a revolving nosepiece with two objective lenses. To switch between them you have to rotate the nosepiece.

## **Changing halogen lamps and fuses (Levenhuk 3ST)**

**CAUTION:** Before changing a lamp or a fuse make sure the microscope is unplugged from a power source. Also, wait until the lamp cools down.

## **Changing the upper illumination source lamp (Levenhuk 3ST)**

Loosen the locking screw on the upper illumination source housing and remove the light socket from within. Replace the old lamp with a new one and put the light socket back. Retighten the locking screw.

## **Changing the lower illumination source lamp (Levenhuk 3ST)**

Loosen the locking screw on the stage and remove the glass. Remove the old lamp from the light socket and replace it with a new one. Put the glass back in its place and retighten the locking screw.

## **Changing the fuse**

Fuse box is located on the bottom of the microscope base. Remove the cover and replace the old fuse with a new one.

## **Care and maintenance**

- After unpacking your microscope and before using it for the first time check for integrity and durability of every component and connection.
- Do not touch the optical surfaces with your fingers. You can use a soft napkin or a cleaning wipe, dipped in ether or absolute alcohol, to clean lenses. We recommend using special optics cleaning fluids from Levenhuk. Do not use any corrosive or acetone-based fluids to clean the optics.
- Abrasive particles, such as sand, should not be wiped off lenses, but instead blown off or brushed away with a soft brush.

- Do not apply excessive pressure when adjusting focus. Do not overtighten the locking screws. Protect the device from sudden impacts and excessive mechanical force.
- Store the device in a dry, cool place away from dust, hazardous acids and other chemicals, away from heaters, open fire and other sources of high temperatures. Do not use the device for lengthy periods of time, or leave it unattended in direct sunlight. Keep the device away from water and high humidity.
- Be careful during your observations, always replace the dust cover after you are finished with observations to protect it from dust and stains.
- If you are not using your microscope for extended periods of time, store the objective lenses and eyepieces separately from the microscope.

## Caution

Never, under any circumstance, look directly at the Sun, another bright source of light or at a laser through this device, as this may cause PERMANENT RETINAL DAMAGE and may lead to BLINDNESS.

Take necessary precautions when using the device with children or people, who have not read and fully understood these instructions.

Do not try to take the device apart on your own. For repairs of any kind, please contact your local specialized service center.

## International Warranty

All Levenhuk telescopes, microscopes, binoculars and other optical products, except for accessories, carry a **lifetime warranty** against defects in materials and workmanship. **Lifetime warranty** is a guarantee on the lifetime of the product on the market. All Levenhuk accessories are warranted to be free of defects in materials and workmanship for **six months** from date of retail purchase. Levenhuk will repair or replace such product or part thereof which, upon inspection by Levenhuk, is found to be defective in materials or workmanship. As a condition to the obligation of Levenhuk to repair or replace such product, the product must be returned to Levenhuk together with proof of purchase satisfactory to Levenhuk.

This warranty does not cover consumable parts, such as batteries.

A Return Authorization (RA) Number must be obtained in advance of return. Contact the local Levenhuk branch to receive the RA number to be displayed on the outside of your shipping container. All returns must be accompanied by a written statement setting forth the name, address and telephone number of the owner, including a description of any claimed defects. Parts or products for which replacement is made will become the property of Levenhuk.

The customer will be responsible for all costs of transportation and insurance to and from Levenhuk or its authorized dealers and will be required to prepay such costs.

Levenhuk will use reasonable efforts to repair or replace any product covered by this warranty within thirty days of receipt. If a repair or replacement will require more than thirty days, Levenhuk will notify the customer accordingly. Levenhuk reserves the right to replace any product that has been discontinued from its product line with a new product of comparable value and function.

This warranty does not apply to any defects or damages resulting from alteration, modification, neglect, misuse, usage of improper power sources, damage in transportation, abuse, or any cause other than normal use, or to malfunction or deterioration due to normal wear.

Levenhuk disclaims all warranties, express or implied, whether of merchantability or fitness for a particular use, except as expressly set forth herein. The sole obligation of Levenhuk under this limited warranty will be to repair or replace the covered product, in accordance with the terms set forth herein. Levenhuk disclaims liability for any loss of profits, loss of information, or for any general, special, direct, indirect or consequential damages which may result from breach of any warranty, or arising out of the use or inability to use any Levenhuk product. Any warranties which are implied and

which cannot be disclaimed will be limited in duration to a term of six months for accessories from the date of retail purchase.

Some states/provinces do not allow the exclusion or limitation of incidental or consequential damages, so the above limitations and exclusions may not apply to you. This warranty gives you specific legal rights, and you may have other rights which vary from state to state or province to province.

Levenhuk reserves the right to modify or discontinue any product without prior notice.

NOTE: This warranty is valid to USA and Canadian customers who have purchased this product from an authorized Levenhuk dealer in the USA or Canada. Warranty outside the USA or Canada is valid only to customers who purchased from an authorized Levenhuk dealer in the specific country or international distributor. Please contact them for any warranty service.

If warranty problems arise, or if you need assistance in using your product, contact the local Levenhuk branch:

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Levenhuk® is a registered trademark of Levenhuk, Inc.
© 2006-2014 Levenhuk, Inc. All rights reserved.
Levenhuk, Inc. 1935 Brandon Court, Suite A-1